

Corning® Erlenmeyer Flasks


Guidelines for Use

CORNING

You can use the fill volumes and speeds shown below as a starting point for optimizing your shaker flask protocol with Corning polycarbonate Erlenmeyer flasks. If you have questions or need assistance, please contact your local Corning office.



	Flask Size	Corning Cat. No.	Recommended Fill Volume (mL)	Recommended Speed (rpm)	Orbital Shaker Size (mm)
CHO-S	125 mL	431143	30	130	25
	250 mL	431144	60	130	25
	500 mL	431145	150	130	25
	1L	431147	350	130	25
	2L	431255	400	130	25
	3L	431252	800	90	25
	5L	431685	2,500	90	25
HEK	125 mL	431143	30	130	15
Sf9	125 mL	431143	30	130	15
	500 mL	431145	200	150	31.75
	1L	431147	300	120 - 150	31.75
	3L	431252	1,000	100 - 150	31.75
	5L	431685	3,500	90 - 100	31.75

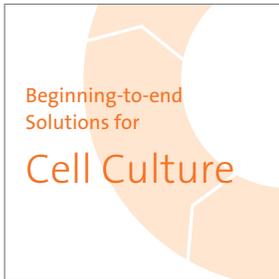

www.corning.com/lifesciences/solutions

At Corning, cells are in our culture. In our continuous efforts to improve efficiencies and develop new tools and technologies for life science researchers, we have scientists working in Corning R&D labs across the globe, doing what you do every day. From seeding starter cultures to expanding cells for assays, our technical experts understand your challenges and your increased need for more reliable cells and cellular material.

It is this expertise, plus a 160-year history of Corning innovation and manufacturing excellence, that puts us in a unique position to offer a beginning-to-end portfolio of high-quality, reliable cell culture consumables.

For more specific information on claims, visit the Certificates page at www.corning.com/lifesciences.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

For additional product or technical information, visit www.corning.com/lifesciences or call 1.800.492.1110. Outside the United States, call +1.978.442.2200 or contact your local Corning sales office.

**Corning Incorporated
Life Sciences**
836 North St.
Building 300, Suite 3401
Tewksbury, MA 01876
t 800.492.1110
t 978.442.2200
f 978.442.2476
www.corning.com/lifesciences

**Worldwide
Support Offices**

ASIA/PACIFIC
Australia/New Zealand
t 61 427286832

China
t 86 21 3338 4338
f 86 21 3338 4300

India
t 91 124 4604000
f 91 124 4604099

Japan
t 81 3-3586 1996
f 81 3-3586 1291

Korea
t 82 2-796-9500
f 82 2-796-9300

Singapore
t 65 6572-9740
f 65 6861-2913

Taiwan
t 886 2-2716-0338
f 886 2-2516-7500

EUROPE

France
t 0800 916 882
f 0800 918 636

Germany
t 0800 101 1153
f 0800 101 2427

The Netherlands
t 020 655 79 28
f 020 659 76 73

United Kingdom
t 0800 376 8660
f 0800 279 1117

**All Other European
Countries**
t 31 (0) 20 659 60 51
f 31 (0) 20 659 76 73

LATIN AMERICA
grupoLA@corning.com

Brasil
t (55-11) 3089-7400

Mexico
t (52-81) 8158-8400

CORNING | **FALCON** | **AXYGEN** | **GOSELIN** | **PYREX**

For a listing of trademarks, visit www.corning.com/clstrademarks.
All other trademarks are the property of their respective owners.