BioMix™ Red

Shipping: On Dry/Blue Ice Catalog numbers

Batch No.: See vial BIO-25006 : 500 x 50 μ l reactions: (10 x 1.25 ml)

Concentration: 2x

BIOLINE A Meridian Life Science® Company Store at -20°C

Storage and stability:

The BioMix Red is shipped on dry/blue ice. On arrival store at -20 °C for optimum stability. Repeated freeze/thaw cycles should be avoided.

Expiry:

When stored under the recommended conditions and handled correctly, full activity of the kit is retained until the expiry date on the outer box label.

Safety precautions:

Please refer to the material safety data sheet for further information.

Routine PCR applications

High throughput

Products suitable for TA cloning

Quality control specifications:

BioMix Red and its components are extensively tested for activity, processivity, efficiency, sensitivity, absence of nuclease contamination and absence of nucleic acid contamination prior to release.

Notes:

Research use only.

BioMix is a Trademark of Bioline Reagents Limited.

Applications

Features

- Convenient pre-mixed, pre-optimized 2x solution
- Premium Tag polymerase suited to a wide range of applications
- Processes fragments up to 5 kb
- Reduced risk of contamination
- Dramatically decreases the time required for reaction set-up
- Reproducible results
- Direct gel loading

Description

BioMixTM Red is a complete ready-to-use 2x reaction mix containing an ultra-stable Tag DNA polymerase. It contains an additional inert red dye that permits easy visualization and direct loading onto a gel. There is no need to add loading buffer as the mix is of sufficiently high density to sink to the bottom of the

BioMix Red has been developed to perform PCR assays of many common genomic and cDNA templates; the user has simply to add water, template and primers. It dramatically reduces the time required to set up reactions, thereby minimizing the risk of contamination. Greater reproducibility is ensured, by reducing the number of pipetting steps that can lead to errors.

BioMix Red has been optimized for a wide variety of templates, however an additional 50mM of MgCl₂ solution is included should any fine adjustments be required.

Components

	500 Reactions
BioMix Red	10 x 1.25 ml
50 mM MgCl ₂ Solution	1.2 ml

Product Citations:

- 1 Ciric, L., et al. FEMS microbial. Let., 353(2): 106-115 (2014).
- Kapralov, M. V., et al. Mol. Boil. Evol. 30(5): 1051-1059 (2013).
- Harrup, L. E., et al. J. Med. Entomol. 49(1): 112-121 (2012).
- Wang, Z., et al. J. Biol. Chem. 286(48): 41359-41367 (2011).
- Price, N. T., et al. J. Biol. Chem. 285(11): 7857-7865 (2010).
- Schultz, J.K., et al. J. Hered. 100(1), 25-33 (2009).
- 7 Kane, N., et al. Endo. J. 150(6), 2882-2888 (2008).
- 8 Rayner, B.S., et al. J. Neurochem. 97(1), 211 (2006).
- Martinez-Cuesta, M.C., et al. Lett. Appl. Microbiol. 40(1), 44 (2005).

BioMix Red Protocol

Reaction Conditions (For a 50 µl reaction)

The optimal conditions will vary from reaction to reaction and are dependent on the system used. Each parameter has to be adjusted individually and some optimization may be required.

BioMix Red	25 µl
Template and Primers	as required
Water (ddH ₂ O)	up to 50 µl

Denature: 94-96 °C

Extension: 70-72 °C Allowing 15-30 seconds per kb

For optimal resolution of PCR products, we recommend the use of Tris-Acetate EDTA (TAE) buffer for gel preparation and electrophoresis.

The Mg²⁺ concentration in the 2x mix is 5 mM (2.5 mM final concentration), this is the optimum concentration for BioMix and should only be adjusted if absolutely necessary. The table below shows the volume of additional MgCl2 to add to a 50 µl reaction to achieve different final concentrations

Final Magnesium concentration required	Volume of 50mM MgCl₂ to add to a 50µl final reaction volume	
2.5 mM	0	
3.0 mM	0.5 μΙ	
3.5 mM	1 µl	

This data is intended for use as a guide only; conditions will vary from reaction to reaction and may need optimization.

Bioline Reagents Ltd UNITED KINGDOM Tel: +44 (0)20 8830 5300

Fax: +44 (0)20 8452 2822

Bioline USA Inc.

Fax: +1 508 880 8993

Bioline GmbH GERMANY Tel: +1 508 880 8990

Tel: +49 (0)337 168 1229 Fax: +49 (0)3371 68 1244 Bioline (Aust) Pty. Ltd AUSTRALIA

Bioline France FRANCE

Meridian Bioscience Asia Pte Ltd SINGAPORE

Tel: +61 (0)2 9209 4180 Fax: +61 (0)2 9209 4763 Tel: +33 (0)1 42 56 04 40 Fax: +49 (0)9 70 06 62 10

Tel: +65 6774 7196 Fax: +65 6774 6441