

FEATURES

Time / calendar display: Hour, minute, seconds, month, date and

Stopwatch: 1 / 100 sec. for 30 min., 1 sec. up to 24 hours. Changeover between 12 - and 24 - hour indications. Note: Depress A to select the desired mode.

NORMAL MODE

(10:09 36)

CALENDAR (è 8)

Hour, minute and seconds are displayed.

Month / Date Keep B depressed. Month, date and day are displayed.

12 - OR 24 - HOUR INDICATION

(a 20:Qi)

(io:09 x)

Select 12 or 24 hour cycle option when setting normal time, hour by advancing figure to "A" / "P" for 12 hour mode or "H" for 24 hour mode

ALARM (<u>1:30 %</u>

Keep C depress to read the alarm time (This illustration shows that the alarm keeps at 7:30AM.)

ALARM ON / OFF

7:30 Alarm engaged.

Reset

Keep B and C depressed at the same time.

STOPWATCH MODE

Depress A to activate this mode B : Start / Stop


C : Split / Split release / Reset

Depress C to reset the stopwatch to "0:0000" before using the stopwatch.

Accumulated elapsed time measurement: B - B - B - B - C



Split time measurement: B - C - C - B - C

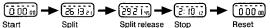


Measurement between two competitors: B - C - B - C - C



Depress A to return to the normal mode.

TIME / CALENDAR SETTING MODE

Depress A three time to activate this mode. Depress C to select the digits to be adjusted (flashing).

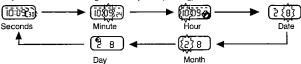

Depress B to reset the seconds to "00" or to set the other digits.

Keep B depressed to advance the digits quickly.

Note: When the seconds count any number from "30" to "59" and C is depressed, one minute is added.

Depress A to return to the normal mode.

ALARM SETTING MODE

(1 } 30 _A)

HOUR

MINUTE

Depress A twice to activate this mode. Depress B to advance the hour.

Depress C to activate this mode. Depress B to advance the minute. Depress A to return to the normal time mode.

HOW TO SET THE HOUR CHIME ON

Depress C, alarm time displayed. Hold It.

Depress A, until all marks above the days of the week appear

The hour chime is on. The stopwatch is set with hour chime.

HOW TO SET THE HOUR CHIME OFF

Depress C, alarm time displayed. Hold it.

Depress A, all marks will disappear, the hour chime is off. No hour Chime.

HOW TO SETTHE SNOOZE FUNCTION ON

When the alarm sounds, depress B, then the alarm will turn off temporarily; Approximately 5 minutes later, the alarm will sound again.

HOW TO SET THE SNOOZE FUNCTION OFF

If you do not need the snooze function, depress C, when the alarm sounds. The alarm will be off, and the snooze function will also be turned off.

If the alarm sounds for more than one minute, without any buttons being depressed, the alarm will automatically turn off and the snooze function will also turn off automatically.